



415 S Airpark Road
Cottonwood, AZ 86326, USA

Phone: (800) 733-0266
Fax: (928) 649-2306



F25/26 Moulton Park Business Centre,
Redhouse Road
Northampton, NN3 6AQ, UK

Phone: 01604 497516
Fax: 01604 497501

DISCOVERY MX10 INCUBATOR



DISCOVERY MX10 PROFESSIONAL INCUBATOR

The Discovery MX10 is designed to incubate up to 8 (10 screw top type) Dipslides, and can be powered from the mains via the adaptor supplied, or via a cigar lighter cable in a vehicle.

Prior to Use:

For mains use, select the applicable plug for your location and clip in the adaptor. The Power input is located on the LH side of the Discovery case. Using the incubator in a vehicle is simply a matter of changing the cable.

A QUICKSET guide point is shown on the temperature control at 37°C and should be used for reference. Each quarter turn of the temperature control equates to 7°C +/- from this point. For example, roughly midway (12.00) would be around 30°C, ideal for L8 dipslide applications.

The actual temperature must always be taken against the thermometer inside the unit, as with all incubators, large variants in ambient temperatures can influence internals slightly. Prior to loading samples, it is recommended that the incubator should be warmed up for a period of 30 minutes to stabilize mains power.

Safety:

The mains power pack features automatic overload protection, and the 12v cigar lighter cable has a 1.5a QB fuse in the large plug end. Please refer any service or repair to Lynd Product Ltd, or your supplier.

Cleaning:

A wipe over with a damp cloth is the best method. Avoid acetone or solvent based cleaners.

Service:

The Incubator is designed for constant long-term operation, and does not contain any user-serviceable parts internally. Please refer to the manufacturer/supplier for repair if required.

Warranty:

This product is under 24-month warranty in normal use. In the event of component failure or defective workmanship, the unit will be repaired or replaced at the company's discretion on an ex-works basis. This comprehensive warranty, which does not affect any statutory rights, applies to the original purchaser only. The date of invoice is deemed to be date of supply. This warranty specifically excludes damage to any part of the Discovery caused by misuse, or any fault caused by incorrect treatment.



End of Life – This product can be Re-manufactured if returned to the supplier/manufacturer – do not dispose of in domestic waste.

Designed & Manufactured in the United Kingdom