



415 S Airpark Road
Cottonwood, AZ 86326, USA

Phone: (800) 733-0266
Fax: (928) 649-2306



F25/26 Moulton Park Business Centre,
Redhouse Road
Northampton, NN3 6AQ, UK

Phone: 01604 497516
Fax: 01604 497501

NUT/NUT

MICROSLIDE® TECHNICAL

DOCUMENT



Index

USE	Page 3
APPLICATION	Page 3
PADDLE AGARS	Page 3
STORAGE/EXPIRATION	Page 3
AGAR VERIFICATION	Page 4
SAMPLING	Page 4
INCUBATION	Page 5
COLONY MEASURING	Page 5
ENUMERATION	Page 5
DISPOSAL	Page 6
IDENTIFICATION	Page 6
GLOSSARY	Page 9

NUT/NUT

CODE: M-NUT/NUT

USE

Isolation and differentiation of Gram (-) enteric bacilli.

APPLICATION

In total coliform testing (TCC), the coliform organisms tested for include: total coliform, fecal coliform, and *E. coli* (*Escherichia coli*). Detection of fecal coliforms (a subset of total coliforms) or *Escherichia coli* (a subset of fecal coliforms) can indicate the potential presence of waterborne pathogens associated with fecal contamination¹.

PADDLE AGARS



Side 1&2: Nutrient-TTC Agar (NUT) – (Color: Yellow) General purpose (relatively non-selective) medium, which will support the growth of a wide variety of organisms. Suitable for cultivation of both aerobes and anaerobes. Aerobic coliform bacteria can be detected by their ability to reduce the TTC dye to a red-colored formozan dye. Bacterial colonies appear as red dots on an otherwise yellow medium.

Note: The Nutrient-TTC agar color is normally light yellow when the agar is cast. After testing, during the incubation phase, the agar may change to a light green color. This color change is a result of either a microbial-induced or chemically-induced pH change in the media. This color change alone does not indicate the presence of microorganisms. Development of red spots or other growth on the agar are an indication of microorganisms.

***Note:** Side 1 of each paddle is marked with an indented laser line.

STORAGE / EXPIRATION

Microslides® should be stored tightly sealed (unopened) in a cool, dry location at room temperature (18 - 25°C; 65 - 77°F). Temperature fluctuations may result in condensation settling at the bottom of the vial, although this does not affect culture properties, it could reduce the shelf-life or cause the agar to separate from the plastic paddle support. Refer to 'Best Before End date' (SEE: BBE stamped on vial).

Avoid sudden temperature changes. Shield from direct sunlight. Do not allow paddles to freeze. Do not store in a refrigerator (~44°F / 10°C) or at temperatures exceeding 80°F; 27°C. Refrigeration may result in water condensation. Discard if paddle agar appears oxidized (darkened from expected color) or if contaminants appear. Expiry applies to medium in its intact container when stored as directed.

AGAR VERIFICATION

¹ United States Pharmacopeial Convention. 2007. The United States pharmacopeia, 31st ed., Amended Chapters 61, 62, 111. The United States Pharmacopeial Convention, Rockville, MD.

These agars have been verified by [EMSL Analytical, Inc.](#) using *E. coli* and *E. faecalis* cultures. Documentation available upon request.

SAMPLING

SURFACE Sampling Protocol

1. Remove the paddle from the vial. Do not touch the agar surfaces.
2. To assure an accurate area recovery, contact the paddle to 20²cm of the surface by contacting the surface twice in separate 10²cm areas.
3. Replace paddle in vial.
4. Incubate.

LIQUID Sampling Protocol


DIRECT IMMERSION PROTOCOL – low viscous liquids

1. Mix liquid test sample.
2. Remove the paddle from the vial. Do not touch the agar surfaces.
3. When taking the sample:
 - a. Pour 40mL of the sample into the vial (to the printed horizontal fill line; see right). Dip the paddle into the 40mL volume liquid in the vial. Maintain a contact time of at least 15 seconds (30 seconds optimal). Both agar surfaces must be completely contacted.
 - b. Or dip the paddle into the sample directly. Maintain a contact time of at least 15 seconds (30 seconds optimal). Both agar surfaces must be completely contacted.
4. Allow excess fluid to drain off both paddle agar surfaces.
5. Replace paddle in vial.
6. Incubate.



SPREAD Protocol – high viscous liquids

1. Mix liquid test sample.
2. Remove paddle from vial. Do not touch the agar surfaces.
3. Holding the contact agar surface on a horizontal plane, deposit volume as a single drop approximately 1cm from the handle boundary (Figure 1).
4. Position a sterile glass rod on the "handle" side of the drop and bring it into contact with the drop creating a meniscus. Drag the glass tube over the paddle agar surface.
5. Replace paddle in vial.
6. Incubate.

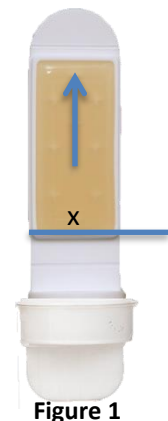


INCUBATION

Incubation of Paddle Growth	Incubation Temperature	Examine at:
Yeast / Mold	25 to 30°C	48 hours up to 120 hours (5 days)
Yeast / Mold	Room Temperature	Up to 7 days
Total Coliform / Bacteria	35 ± 2°C	24 to 48 hours
Total Coliform / Bacteria	Room Temperature	Up to 5 days

Note: Incubation of bacteria after 48 hours may produce confluent growth making enumeration more difficult.


COLONY MEASURING

Each Microslide® paddle has molded media attachment points that are 4mm in length (point-to-point). This feature provides a useful guidepost to estimating nearby colony size.

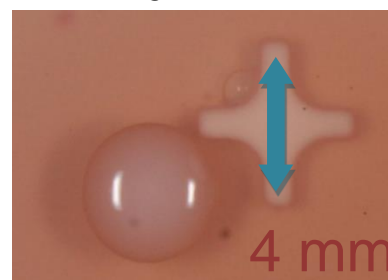

ENUMERATION

Bacteria CFU/mL

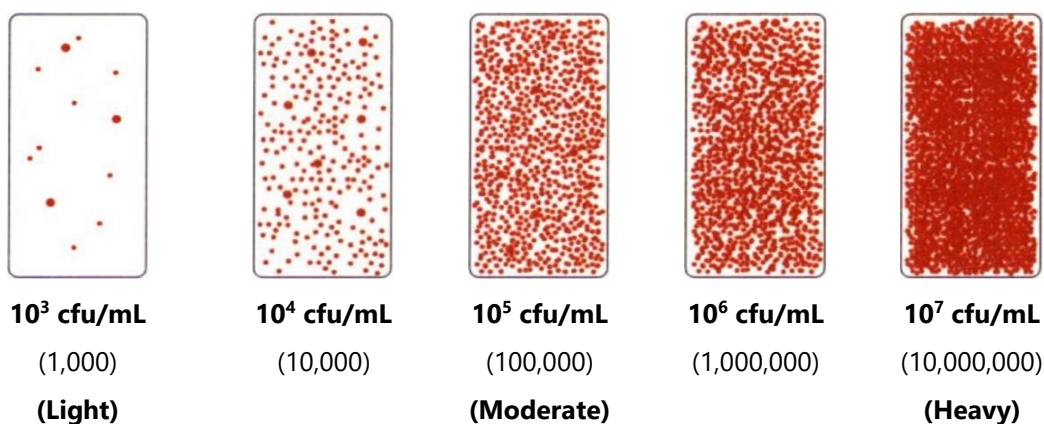


Note: Estimation of lower counts is possible, but statistically difficult to justify. Use Light, Moderate and Heavy for Mold growth and surface testing.

DISPOSAL

Make a 1:9 dilution of household bleach (5.25% sodium hypochlorite solution). Twist and remove Microslide® paddle from vial. Fill vial with 40mL diluted hypochlorite solution (to fill-line). Allow 15-minute contact time. Discard bleach solution. Replace paddle in vial and dispose. Alternatively, loosen cap and microwave for 30 seconds, autoclave, or incinerate.

IDENTIFICATION

Organism	Nutrient-TTC (NUT)
<i>Aspergillus niger</i>	 <p>Growth: +++ Colony: Granular, jet black conidia with yellow/gray hyphae, 3-5++cm</p>
<i>Bacillus spp.</i>	 <p>Growth: +++ Colony: Opaque with dark center (bullseye), irregular, raised, lobate (wrinkled), 2-4mm+</p>
<i>Candida albicans</i>	 <p>Growth: +++ Colony: Cream, CVEG, 1-2mm</p>
<i>E. coli</i>	 <p>Growth: +++</p>

Enterobacter aerogenes

Colony: Yellow/Orange/Red, CVEG, 2-4mm




Growth: +++

Colony: Maroon/red with transparent margin, CVEG, 0.1-0.5mm

Enterococcus spp.

INHIBITED

Klebsiella spp.

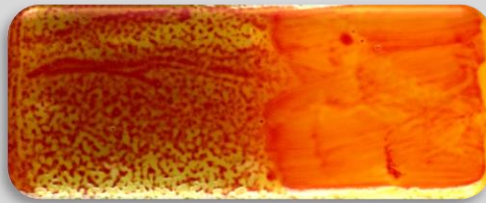

Growth: +++

Colony: Amber/Red, spreading, 0.5-1.0mm

Proteus spp.



Growth: +++

Colony: Maroon/red with dark red center and transparent margin, irregular, glistening (swarming-transparent field), raised, undulate, 1-4mm

Pseudomonas aeruginosa



Growth: +++

Colony: Maroon/red with transparent margin, circular to irregular, raised, entire, 1-2mm

Pseudomonas fluorescens



Growth: +++

Salmonella enteritidis


Colony: Clear/colorless with grey/dark center, translucent edges, irregular/spreading to confluent, 2-4mm



Growth: +++

Colony: Red, FED, 0.5-1.0mm


Serratia spp.

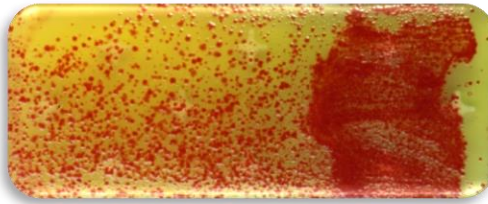

Growth: ++

Colony: Red, FED, 0.5-1.0mm

Shigella spp.

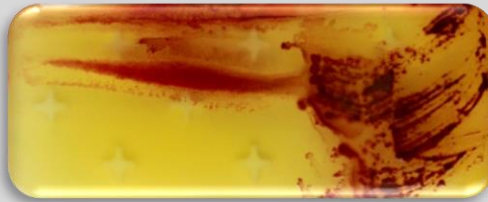

Growth: +

Colony: Maroon/red, CVEG, 0.5-1.0mm

Staphylococcus aureus



Growth: +

Colony: Red, FED, 0.5-1.0mm

Streptococcus spp.



Growth: ++

Colony: Maroon/red, CVEG, 0.1-0.5mm

Streptomyces griseus



Growth: +
Colony: Yellow, FED, 0.5-1.0mm

Gram (+) Bacteria

PARTIAL TO COMPLETE INHIBITION

GLOSSARY

CVEG..... Convex, Entire, Glossy

FED..... Full, Entire, Dull

Gram..... Gram reaction